


Carousel Stand

取扱説明書

この度は本製品をお買い上げいただき、誠にありがとうございました。
ご使用前に本書をよくお読みの上、正しくお使いください。

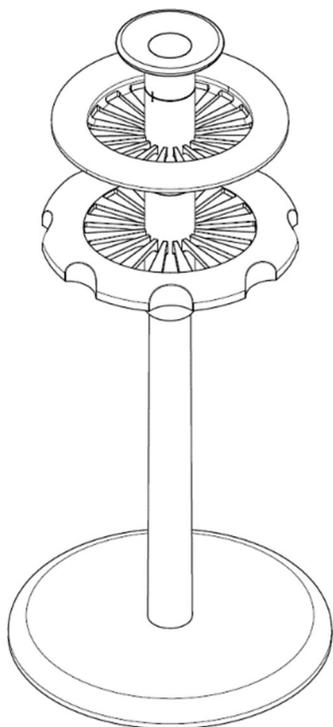

適合機種

Nichipet Air
Nichipet Premium
Nichipet Premium LT
NICHIRYO Le

付属品

ご使用前に下記の付属品がすべて揃っているかご確認ください。


品名	数量
アームホルダー	1
ボディホルダー	1
シャフト	1
ハンドル	1
支柱	1
台座	1
取扱説明書(本紙)	1

使用上の注意事項



- ✓ 本製品はピペットを保管する以外の用途では使用しないでください。
- ✓ ピペットはボディホルダーの凹部に合わせて掛けて下さい。
- ✓ 本製品はピペット8本まで使用可能です。それ以上の本数は使用しないでください。ピペットの落下、スタンド倒れの原因となります。

組立方法

- ①支柱を台座に差し込み、時計方向に回転させ、止まるまで締め付けます。
- ②アームホルダーとボディホルダーを組み合わせた状態で、シャフトを差し込み時計方向に回転させ、止まるまで締め付けます。【図1】
- ③ハンドルとアームホルダーのマークを合わせ、上からハンドルを両サイドの固定穴が見えなくなるまで強く押し込みます。【図2】

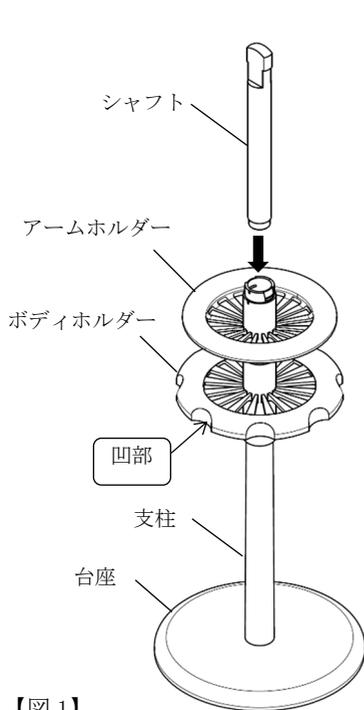

【図1】

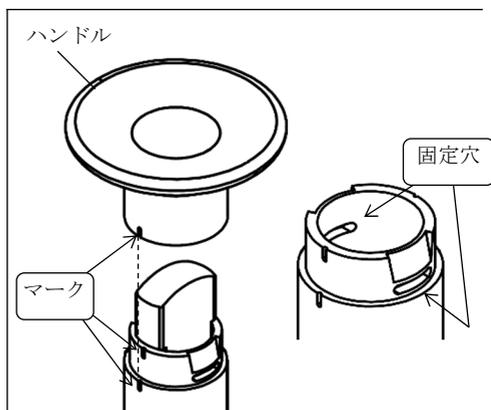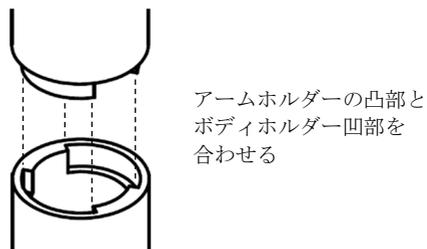

【図2】

販売元

株式会社 **ニチリョー**

Webサイト

<https://www.nichiryo.co.jp/>

Eメール

info@nichiryo.co.jp


●お問い合わせはWebサイトの「お問い合わせフォーム」またはEメールでお願い致します。

SB-NARSTD002EJ

Carousel Stand

User's Manual

Thank you very much for purchasing carousel stand.
Please read this manual carefully before using this product.

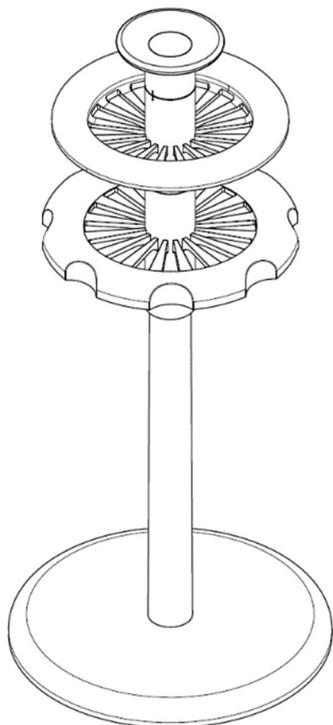

Compatibility

Nichipet Air
Nichipet Premium
Nichipet Premium LT
NICHIRYO Le

Standard component

When unboxing the package, check that all of the following items are included.


Item	QTY
Finger rest Holder	1
Body holder	1
Shaft	1
Handle	1
Column	1
Base	1
User's Manual	1

Safety Precautions


- ✓ Do not use this product for any purpose other than storing the pipette.
- ✓ Hang the pipette in the concave part of the body holder.
- ✓ This product can be used with up to 8 pipettes. Please do not use for more than 8 pipettes. This may cause the pipette to drop or the stand to fall down.

Assembly method


- ① Insert the column into the base, rotate it clockwise, and tighten until it stops moving.
- ② Insert the shaft into the finger rest holder and the body holder combined. Turn clockwise and tighten until it stops moving. (Fig.1)
- ③ Align the mark on the handle and finger rest holder, and fix the handle from above on both sides. Press the handle firmly until the fixing holes to be covered thoroughly. (Fig.2)

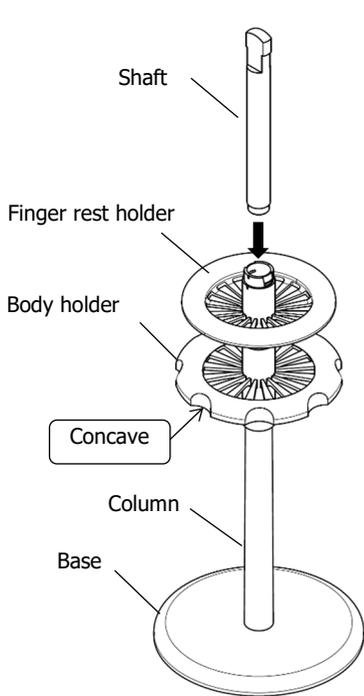

【Fig.1】

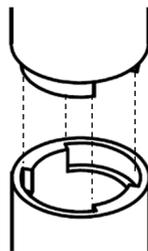

Combine the convex part of the finger rest holder and the concave part of body holder.

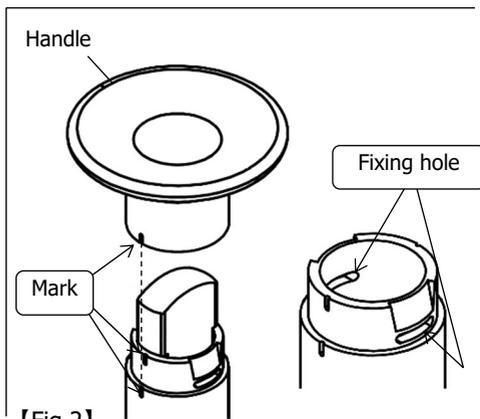

【Fig.2】

Manufacturer :

NICHIRYO CO.,LTD.



Head office & Factory

2760-1 Nishikata, koshigaya-Shi, Saitama 343-0822, Japan
 TEL: +81-48-989-1301 FAX: +81-3-6273-7944

Tokyo office (Overseas Sales Department)

PMO-Kandaiwamoto-cho II 4F,2-2-4 Iwamoto-cho, Chiyoda-ku,
 Tokyo 101-0032, Japan
 TEL: +81-3-5829-4539 FAX: +81-3-5829-4504

URL: <https://www.nichiryo.co.jp/en/>
 E-mail: i-sales@nichiryo.co.jp

SB-NARSTD002EJ