


Multi Rotary Stand

取扱説明書

この度は本製品をお買い上げいただき、誠にありがとうございました。
ご使用前に本書をよくお読みの上、正しくお使いください。

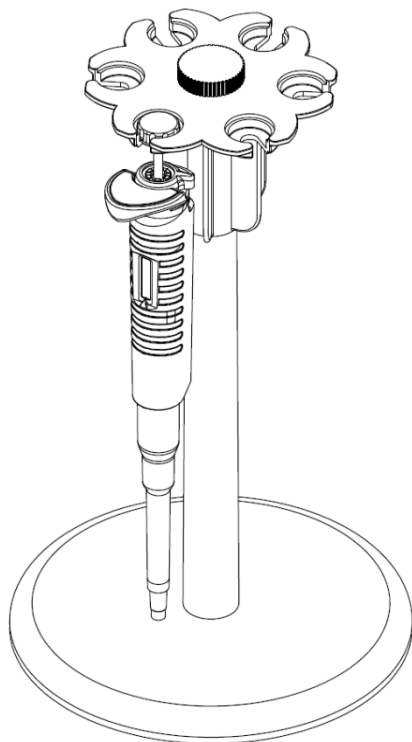

適合機種

Nichipet EX III
Nichipet Premium
Nichipet Premium LT
Nichipet EX II
Nichipet EX II MULTI
Nichipet EX Plus II

付属品

ご使用前に下記の付属品がすべて揃っているかご確認ください。

品 名	数 量
シャフト	1
トッププレート	1
支柱	1
台座	1
取扱説明書(本紙)	1

使用上の注意事項

- ✓ 本製品はピペットを保管する以外の用途では使用しないでください。
- ✓ ピペットはトッププレートの凹部にプッシュボタンを掛けてご使用下さい。



組立方法

- ① 支柱を台座に差し込み、時計方向に回転させ、止まるまで締め付けます。

【図 1】

- ② トッププレートとシャフトを組み合わせた状態で、支柱に差し込みシャフトを時計方向に回転させ、止まるまで締め付けます。

【図 2】

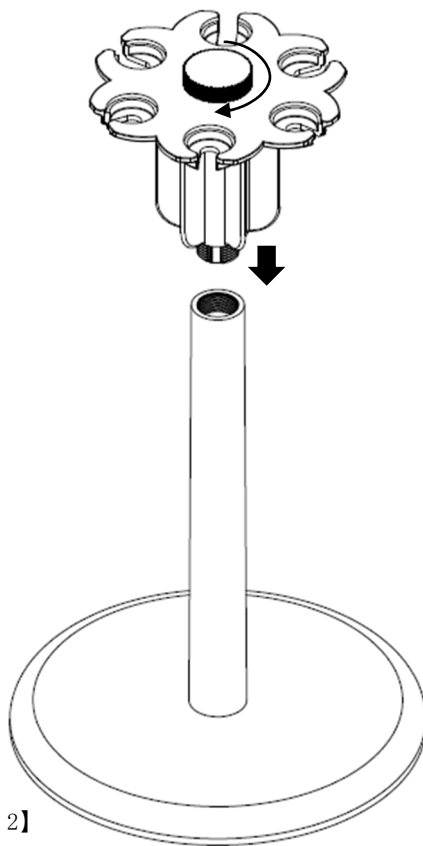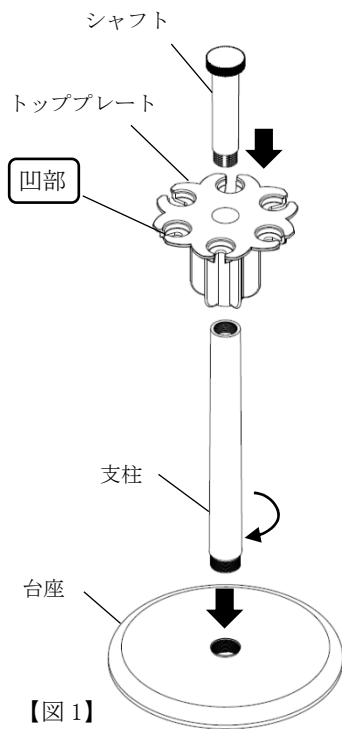

販売元

株式会社 **ニチリョー**

Webサイト <https://www.nichiryo.co.jp/>

Eメール info@nichiryo.co.jp


●お問い合わせはWebサイトの「お問い合わせフォーム」またはEメールでお願い致します。

SB-MLTSTD001EJ

Multi Rotary Stand

User's Manual

Thank you very much for purchasing Multi Rotary Stand.
Please read this manual carefully before using this product.

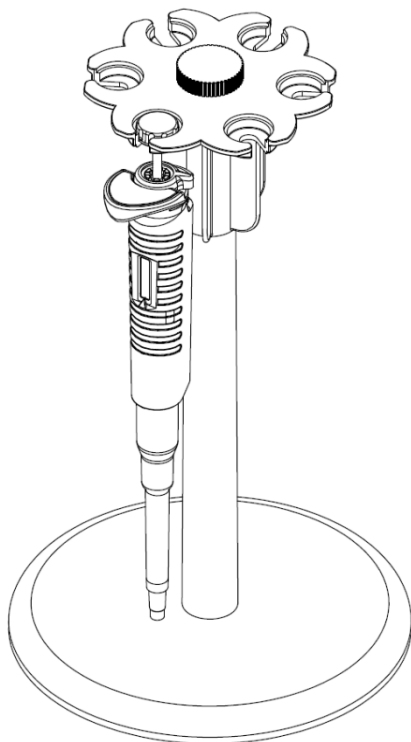

Compatibility

Nichipet EX III
Nichipet Premium
Nichipet Premium LT
Nichipet EX II
Nichipet EX II MULTI
Nichipet EX Plus II

Standard component

When unboxing the package, check that all of the following items are included.

Item	QTY
Shaft	1
Top plate	1
Column	1
Base	1
User's Manual	1

Safety Precautions

- ✓ Do not use this product for any purpose other than storing the pipette.
- ✓ Hang the push button of the pipette in the concave part of the Top Plate.


Assembly method

- ① Insert the Column into the Base, rotate it clockwise, and tighten until it stops moving.


【Figure 1】

- ② Insert the combined parts that Shaft & Top plate into the column. Rotate the shaft clockwise and tighten until it stops moving.

【Figure 2】

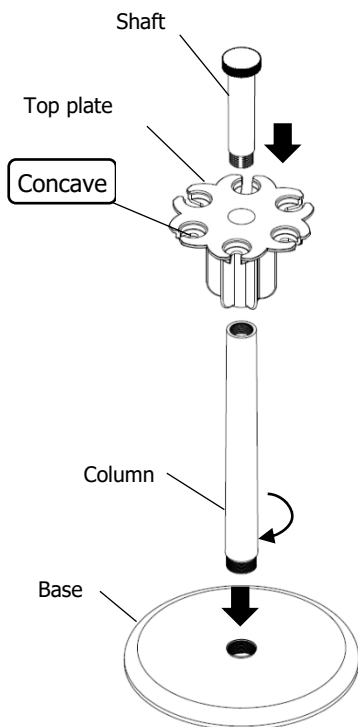

【Figure 1】

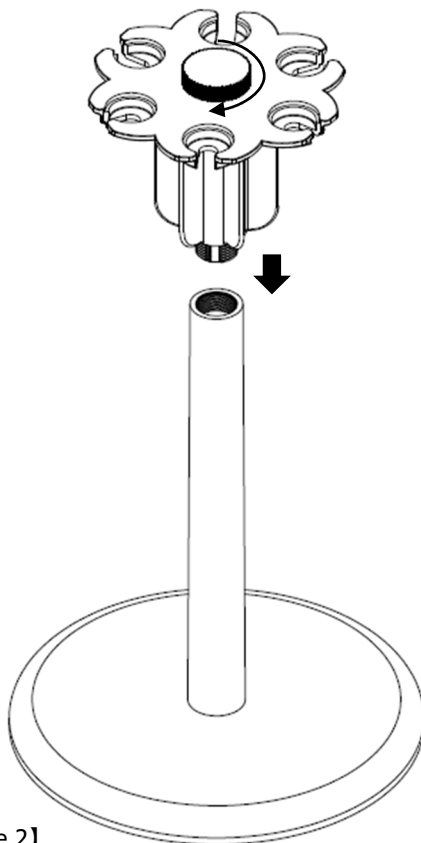

【Figure 2】

Manufacturer :

NICHIRYO CO.,LTD.



Head office & Factory

2760-1 Nishikata, koshigaya-Shi, Saitama 343-0822, Japan

TEL: +81-48-989-1301 FAX: +81-3-6273-7944

Tokyo office (Overseas Sales Department)

PMO-Kandaiwamoto-cho II 4F,2-2-4 Iwamoto-cho, Chiyoda-ku,

Tokyo 101-0032, Japan

TEL: +81-3-5829-4539 FAX: +81-3-5829-4504

URL: <https://www.nichiryo.co.jp/en/>

E-mail: i-sales@nichiryo.co.jp

SB-MLTSTD001EJ