


Standard component

- Silicone tube set ···1 pc
- Filling funnel ···1 pc
- Nozzle cap ···1 pc

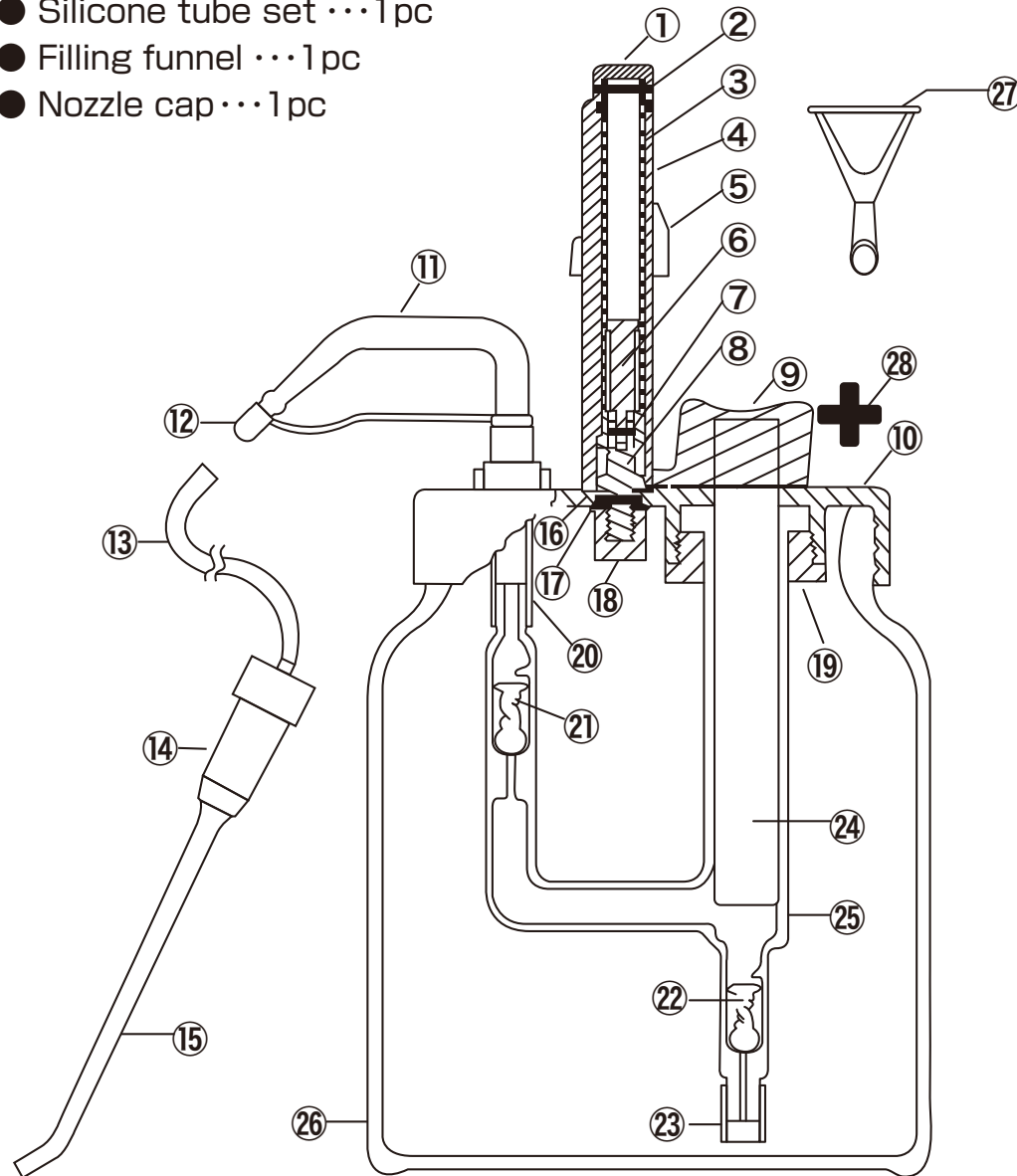

- [ 1 ] Post knob (PP)
- [ 2 ] Post knob screw (Chrome plated brass)
- [ 3 ] Post holder assembly (Chrome plated brass)
- [ 4 ] Post with scale (PC)
- [ 5 ] Volume indicator with screw (PEEK)
- [ 6 ] Operating rod (Chrome plated brass)
- [ 7 ] Spring pin (SUS304)
- [ 8 ] Screw (Chrome plated brass)
- [ 9 ] Plunger assembly (Pyrex, Glass & PP)
- [ 10 ] Bottle cap (PPS)
- [ 11 ] Dispense nozzle (Pyrex glass)
- [ 12 ] Dispense nozzle cap (PP)
- [ 13 ] Silicon tube
- [ 14 ] Tip holder (PP)
- [ 15 ] Glass tip (Glass)
- [ 16 ] Spring pin (SUS304)
- [ 17 ] Washer (PTFE)
- [ 18 ] Lock nut for post holder (PTFE)
- [ 19 ] Lock nut for barrel assembly
- [ 20 ] Barrel/nozzle joint tube (PTFE)
- [ 21 ] Upper screw valve (Glass)
- [ 22 ] bottom screw valve (Glass)
- [ 23 ] Inlet tube (PTFE)
- [ 24 ] Plunger (Glass)
- [ 25 ] Barrel assembly (Pyrex)
- [ 26 ] Amber bottle (Borosilicate glass)
- [ 27 ] Filling funnel (PP)
- [ 28 ] Cap