


Volume adjustment flow

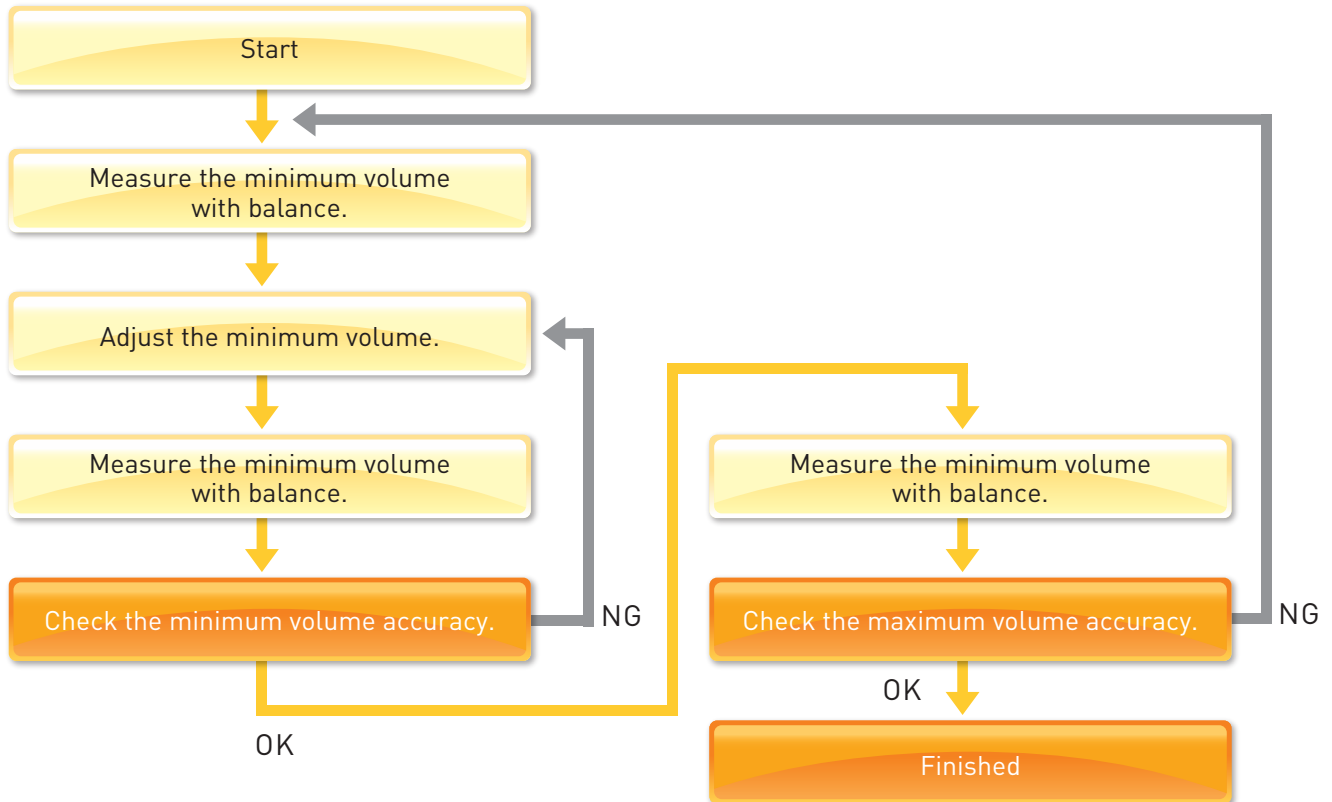

*The minimum volume : The minimum setting volume in each volume range.

*The maximum volume : The maximum setting volume in each volume range.

Reference for volume adjustment by the push button scale (ul)

Type	Correction value for 1/2 scale	Correction value for one scale	Correction value for five scales	Correction value for ten scales
NPLO-2	0.0013	0.0026	0.013	0.026
NPLO-10	0.0065	0.013	0.065	0.132
NPLO-20	0.013	0.026	0.131	0.263
NPLO-100	0.067	0.133	0.665	1.330
NPLO-200	0.132	0.264	1.322	2.644
NPLO-1000	0.659	1.329	6.597	13.194
NPLO-5000	3.332	6.663	33.316	66.633
NPLO-10ML	6.696	13.393	66.966	133.932

Note: Above values are theoretical values. Accuracy and precision should be calculated by actual measured volumes.