


Volume adjustment flow

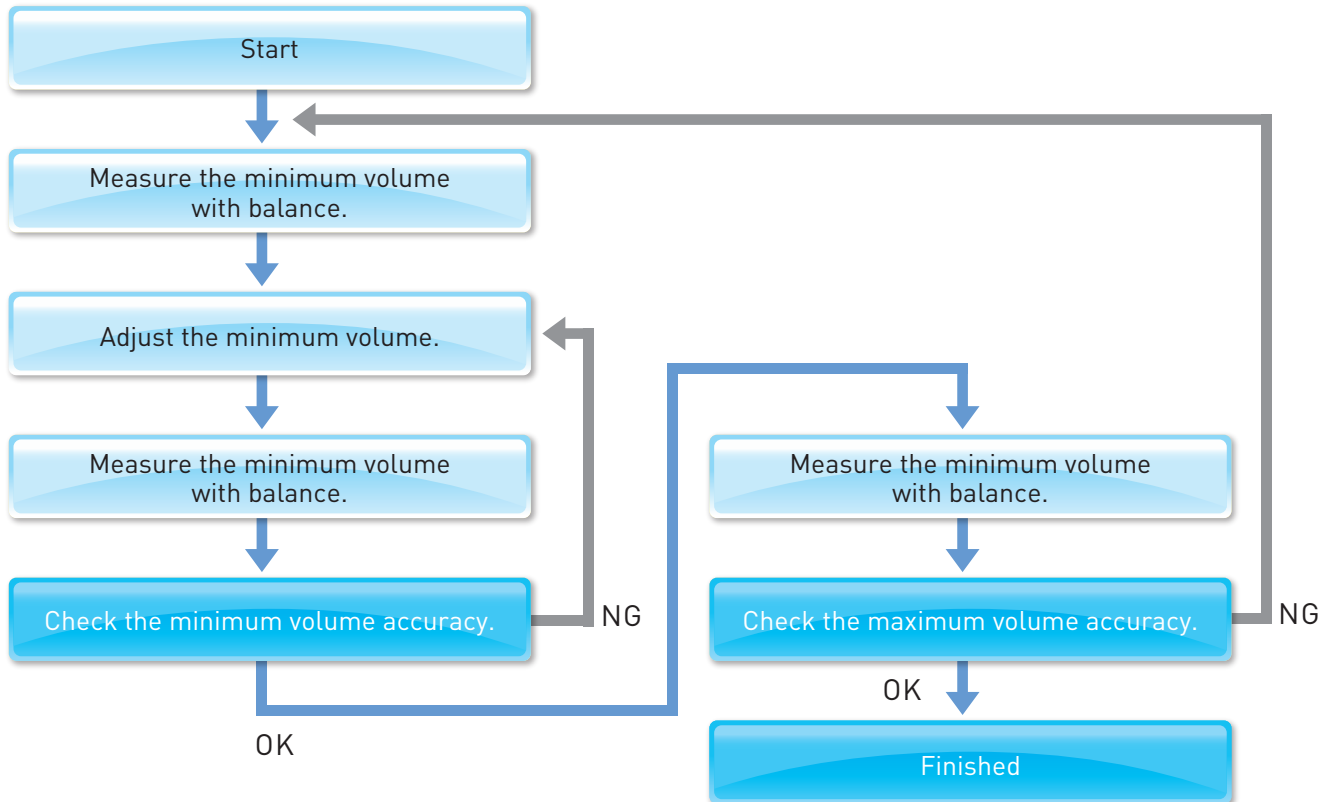

*The minimum volume : The minimum setting volume in each volume range.

*The maximum volume : The maximum setting volume in each volume range.

Reference for volume adjustment by the push button scale (ul)

Type	Correction value for one scale	Correction value for two scale	Correction value for five scales	Correction value for ten scales
NPP-2	0.0015	0.0029	0.0073	0.0145
NPP-10	0.0077	0.0154	0.0385	0.0769
NPP-20	0.012	0.025	0.062	0.125
NPP-100	0.078	0.156	0.39	0.78
NPP-200	0.13	0.25	0.63	1.26
NPP-1000	0.79	1.58	3.96	7.91
NPP-5000	8	16	40	79.9
NPP-10000	8	16.1	40.2	80.3

Note: Above values are theoretical values. Accuracy and precision should be calculated by actual measured volumes.